

How to Add a New Label Printer to StickerBot

by Shannon Patridge | Feb 23, 2022 | Setup | 0 comments



Add a new Dymo, Zebra, or Brother label printer to your Toxilab setup.

Install the new label printer

Find, download, and install the driver for the [Dymo](#), [Zebra](#), or [Brother](#) label printer. It may come on a disc or a download page online. Toxilab and StickerBot do not support any other printers.

Set the new printer as default

Find your network printers via Settings/Bluetooth & devices/Printers & scanners.

Find "Let Windows manage my default printer" and **set this to OFF**.

Locate the printer on the list by name/model and click > to open properties. **Click Set as default**.

Select and test

Open StickerBot and select the channel for your printer.

Send a test job. Troubleshoot if no test label prints.

Search

Recent Posts

Post 2

Hello world!

Recent Comments

Legal

- [Privacy Policy](#)
- [Terms & Conditions](#)
- [Toxilab Community](#)

Contact Us

e: hello@toxilab.com

p: [855.723.7626](tel:855.723.7626)

Toxilab Operations
P.O. Box 931
Getzville, NY 14068

Get Connected

